RotoPol-11

RotoPol-15

Manual No.: 15267001

Date of Release: 01.09.1997

Instruction Manual

RotoPol-11/-15
Instruction Manual

Table of Contents	Page
User's Guide	1
Reference Guide.....	13
Quick Reference Guide.....	22

Always state *Serial No* and *Voltage/frequency* if you have technical questions or when ordering spare parts. You will find the Serial No. and Voltage on the type plate of the machine itself. We may also need the *Date* and *Article No* of the manual. This information is found on the front cover.

The following restrictions should be observed, as violation of the restrictions may cause cancellation of Struers legal obligations:

Instruction Manuals: Struers Instruction Manual may only be used in connection with Struers equipment covered by the Instruction Manual.

Service Manuals: Struers Service Manual may only be used by a trained technician authorised by Struers. The Service Manual may only be used in connection with Struers equipment covered by the Service Manual.

Struers assumes no responsibility for errors in the manual text/illustrations. The information in this manual is subject to changes without notice. The manual may mention accessories or parts not included in the present version of the equipment.

The contents of this manual is the property of Struers. Reproduction of any part of this manual without the written permission of Struers is not allowed.

All rights reserved. © Struers 1997.

Struers A/S
Valhøjs Allé 176
DK-2610 Rødovre/Copenhagen
Denmark
Telephone +45 36 70 35 00
Fax +45 38 27 27 01

RotoPol-11/-15

Safety Precaution Sheet

To be read carefully before use

1. The operator(s) should be fully instructed in the use of the machine, any connected equipment and accessories and the applied consumables according to the relevant Instruction Manuals.
2. The machine must be placed on a safe and stable table with an adequate working height. All functions on the machine and any connected equipment must be in working order.
3. Operators should ensure that the actual voltage corresponds to the voltage on the back of the machine. The machine must be earthed. Follow the local regulations.
4. Connect only to cold water. Make sure that the water connections are leak-proof and that the water outlet is working.
5. The connection to the mains water supply should be turned off if the machine is left unattended.
6. Establish efficient exhaust from the working area if consumables containing alcohol, e.g. blue DP-Lubricant, are being used.
7. Make sure that the grinding or polishing discs are correctly mounted, the driving pins engaged with the holes in the turntable.
8. Keep clear of the rotating disc and the specimen mover during operation. While grinding manually, be careful not to touch the grinding paper. Do not drop the specimen while grinding.
9. Never try to move the specimen mover into upright position while specimens are prepared.
10. If you observe malfunctions or hear unusual noises - stop the machine and call technical service.

The equipment is designed for use with consumables supplied by Struers. If subjected to misuse, improper installation, alteration, neglect, accident or improper repair, Struers will accept no responsibility for damage(s) to the user or the equipment.

Dismantling of any part of the equipment, during service or repair, should always be performed by a qualified technician (electromechanical, electronic, mechanical, pneumatic, etc.).

RotoPol-11/-15
Instruction Manual

User's Guide

Table of Contents	Page
1. Getting Started	
Checking the Contents of Packing.....	2
RotoPol-11 (ROT11).....	2
RotoPol-15 (ROT15).....	2
Recirculation Cooling Unit, Optional (TRECI).....	2
Recirculation Cooling Unit Connector, Optional (ROPRE).....	2
Getting Acquainted with RotoPol-11/-15.....	3
Back of RotoPol.....	3
Unpacking RotoPol.....	4
Placing RotoPol.....	4
Supplying Power.....	4
Supplying Water.....	4
Connecting the Water Inlet.....	4
Connecting the Water Outlet.....	5
Software Settings.....	6
Setting the Language.....	6
2. Basic Operation	
Using the Controls.....	7
The Front Panel of RotoPol-11.....	7
The Front Panel of RotoPol-15.....	7
Front Panel Controls.....	8
General Use of Keys.....	9
Acoustic Signals.....	9
Reading the Display.....	9
Routine Operations.....	10
Mounting a Disc.....	10
Starting the Disc.....	10
Stopping the Disc.....	10
Selecting Water and Speed.....	10
Settings with a RotoCom Memory Unit.....	11

1. Getting Started

Checking the Contents of Packing

RotoPol-11 (ROT11)
RotoPol-15 (ROT15)

In the packing box you should find the following parts:

- 1 RotoPol-11/-15
- 1 Mains cable
- 1 Splash ring
- 1 Inlet hose 3/4" x 2 m
- 1 Filter gasket
- 1 Gasket $\varnothing 11/r24$ x 1.5
- 1 Reduction ring with gasket
- 1 Outlet hose r32 x 3 m
- 2 Hose clamps
- 1 Allen key 5 mm
- 1 Set of Instruction Manuals

Recirculation Cooling Unit,
Optional (TRECI)

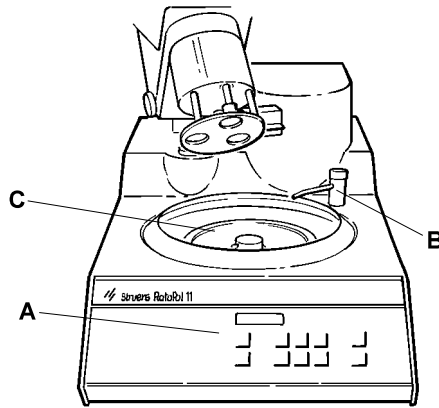
- 1 TRECI complete
- 1 Outlet angle
- 1 Funnel
- 2 Disposable plastic containers (TREPO)

Recirculation Cooling Unit
Connector, Optional (ROPRE)

- 1 Connection box
- 1 Hexagon nipple
- 1 Non-return valve with elbow and hexagon nipple
- 1 Reduction piece, large
- 1 Reduction piece, small
- 1 Hose clamp
- 1 Additive for Cooling Fluid, 1 l (ADDUN)
- 1 Diagram

Getting Acquainted with RotoPol-11/-15


Take a moment to familiarise yourself with the location and names of all the RotoPol components:

RotoPol with specimen mover RotoForce-1 (optional)

- A** Front panel control(s)
- B** Water tap with sprinkler tube
- C** Turntable

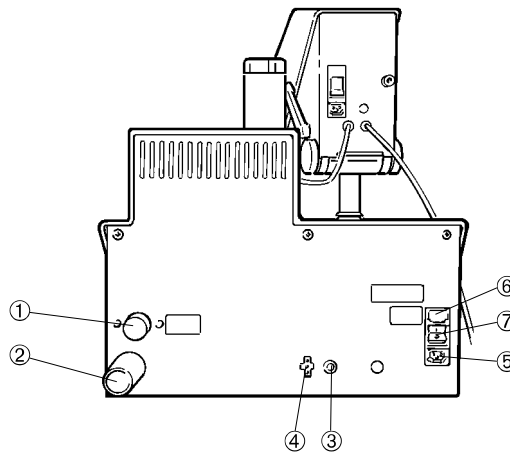Back of RotoPol

- ① Water inlet tube
- ② Water outlet tube
- ③ Socket for specimen mover connection
- ④ Socket for recirculation cooling unit connection
- ⑤ Mains connection
- ⑥ Fuse
- ⑦ Main switch

Unpacking RotoPol

Detach RotoPol from the bottom of the packing case by removing the four screws from below.

Placing RotoPol

Place RotoPol on a steady table with an adequate working height. The machine must be close to the power supply, water mains and water outlet facilities. If water recirculation is used, there must be room under the table for the Recirculation Cooling Unit (TRECI).

Supplying Power

The mains cable of RotoPol is either equipped with a standard electric male plug, type “F” or with a male flat blade attachment plug. If the type of plug delivered with the equipment has not been approved for local use, the plug must be replaced with an approved type. The leads have to be connected as following:

Yellow/green: earth
Brown: phase
Blue or Black: neutral

IMPORTANT

Check that the mains voltage corresponds to the voltage stated on the type plate on the back of the machine.

- Insert the female plug of the power cable into the socket on the back of the machine.
- Connect the male plug to a grounded electric outlet.

DANGER!

The machine must be earthed

Supplying Water

Water for wet grinding may be supplied from the water mains or a Recirculation Cooling Unit.

Connecting the Water Inlet

- Mount the straight end of the pressure hose onto the water inlet tube (upper) on the back of RotoPol (see drawing on the previous page):
 - Insert the filter gasket in the coupling nut with the flat side against the pressure hose.
 - Tighten the coupling nut completely.

- Mount the other end of the pressure hose on the water mains tap for cold water:
 - Mount the reduction ring with gasket on the water mains tap, if necessary.
 - Introduce the gasket and tighten the coupling nut completely.

IMPORTANT
Connect only to cold water.

Connecting the Water Outlet

- Locate the end of the drain hose where a short length of the steel reinforcement has been stripped away. Mount the stripped end of the hose onto the water outlet tube. Use the included hose clamp for fastening.
- Lead the other end of the drain hose to the water outlet. Arrange the tube so that it slopes throughout its length. Shorten the hose, if necessary. The drain hose should be attached to the water outlet system if possible.

REMEMBER
Avoid sharp bends in the drain hose and make sure that the hose slopes throughout its length.

Software Settings

- Switch on the power at the main switch. The following display will appear briefly:

- ROTOPOL-11 -
VERSION X.Y

- ROTOPOL-15 -
VERSION X.Y

- The operation menu now appears:

WATER OFF 150RPM

Setting the Language

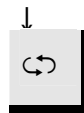
WATER OFF 150RPM

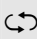
Example screen..


The MENU  key activates the set-up menu.


LANGUAGE
ENGLISH


Press SELECT  to choose between English, German, French or Japanese

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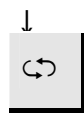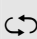Press  until you reach the language you prefer.

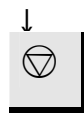LANGUE
FRANCAIS

Press STOP  to accept the choice.

The Operation Menu will appear in the language you have chosen.

EAU OFF 150 T/M

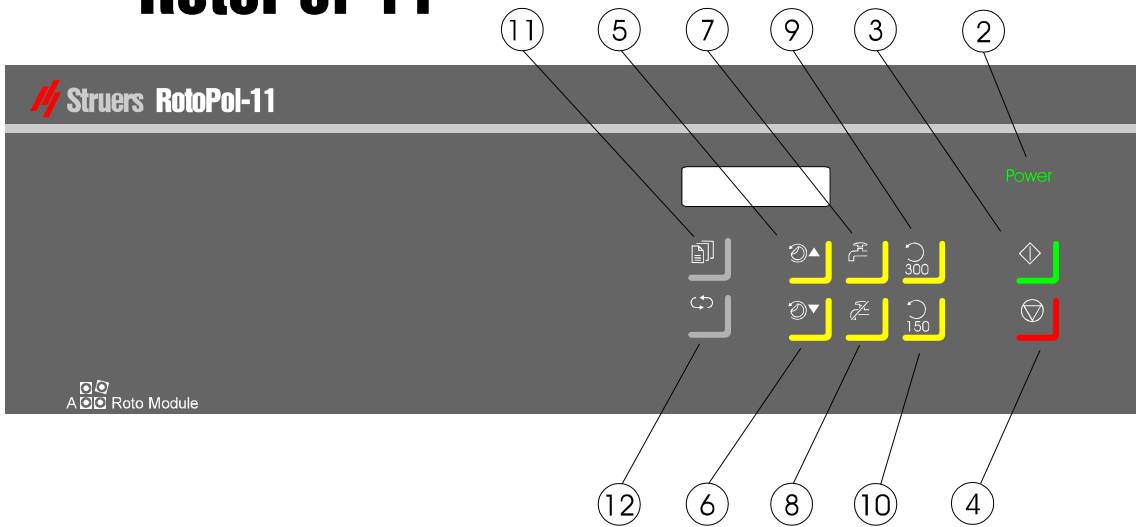
Remember to set the same language on all connected RotoSystem modules. Setting the Communication: see Advanced Operation Section.

2. Basic Operation

Using the Controls

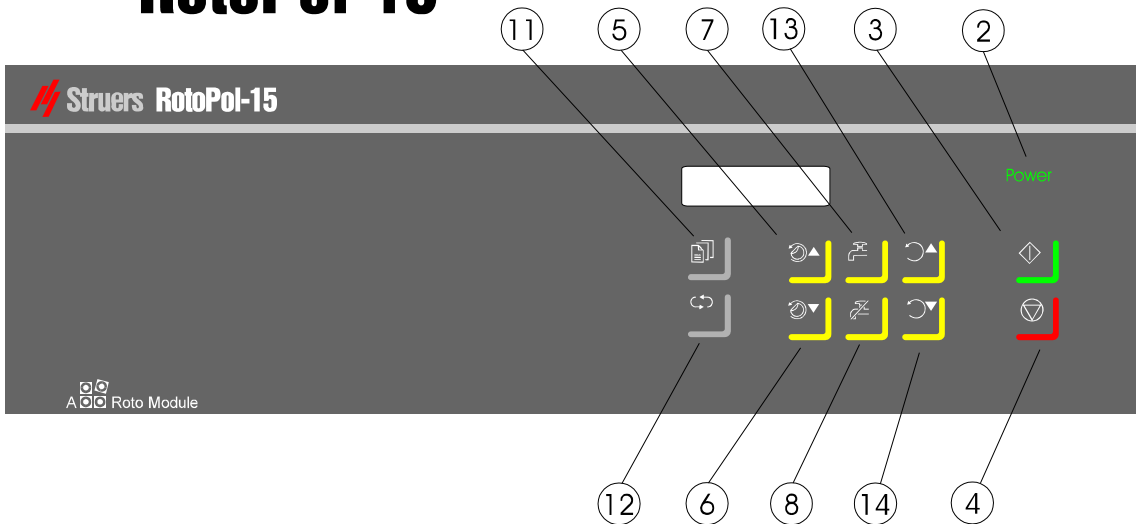
The Front Panel of RotoPol-11

RotoPol-11

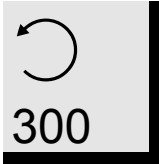
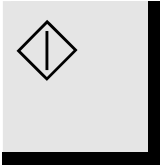
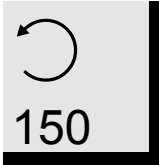
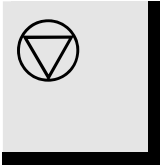
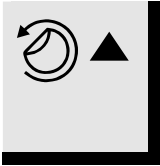
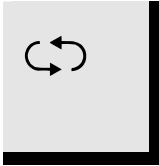
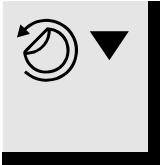
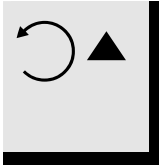
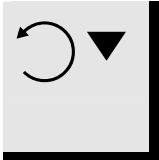
The Front Panel of RotoPol-15

RotoPol-15

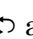
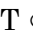
Front Panel Controls

Communication is set to STANDARD mode

Name	Key	Function	Name	Key	Function
① MAIN SWITCH		The main switch is located on the back of the cabinet.	⑧ WATER OFF		Select WATER OFF.
② POWER INDICATOR		Lights up when the main power is on.	⑨ SPEED 300 RPM		Select 300 RPM. (RotoPol-11)
③ START		Starts the disc and any connected RotoSystem equipment set in Auto mode.	⑩ SPEED 150 RPM		Select 150 RPM. (RotoPol-11)
④ STOP		Stops the disc and any connected RotoSystem equipment set in Auto mode. Exits the set-up menu.	⑪ MENU		Activates the set-up menu or moves down in the menu structure.
⑤ DISC		Scrolls up in the operation menu.	⑫ SELECT		Changes values in the set-up menu.
⑥ DISC		Scrolls down in the operation menu.	⑬ SPEED		Increase the speed (RotoPol-15)
⑦ WATER ON		Select WATER ON: press the key briefly. Manual Dosing: keep pressing the key. Release the key to stop the water.	⑭ SPEED		Decrease the speed (RotoPol-15)

General Use of Keys

- WATER ON , WATER OFF  and SPEED  can only be selected freely at User Select.
- MENU , SELECT  and DISC  can only be activated when the machine has been stopped.
- SELECT  can only be activated when MENU  is selected.

Acoustic Signals

- When pressing a key, a short beep indicates that the command has been accepted, whereas a long beep indicates that the key cannot be activated at the moment.

Reading the Display

WATER ON 300RPM
GRINDING

- The first line on the screen shows the setting of parameters WATER and SPEED.
- The second line, if any, shows a selected pre-set mode (GRINDING or POLISHING).

Routine Operations

Mounting a Disc

Place the disc on the turntable and turn it until the three driving pins engage with the holes in the turntable.

Starting the Disc

Press START . The disc starts rotating. The selected functions are executed, and connected RotoSystem modules start operating.

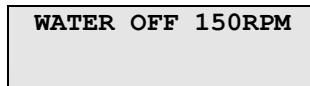
Stopping the Disc

Press STOP . The disc stops rotating. The selected functions and connected RotoSystem modules stop.

Selecting Water and Speed

You can either select water and speed yourself or press DISC   to select one of the pre-set GRINDING and POLISHING modes.

Setting to User Select

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
The User Select screen has an empty line at the bottom.

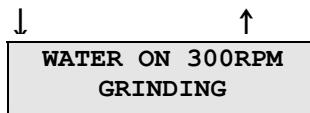
- Press  or  to select WATER ON or WATER OFF.
- Press SPEED  150 or  300 to select the speed. (RotoPol-11)
- Press SPEED  or  to adjust the speed. (RotoPol-15)

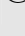
Press DISC  or  to change between the settings.

Setting to GRINDING

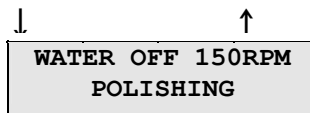
The standard setting GRINDING is pre-set to WATER ON and 300RPM.

Press DISC  or  to change between the settings.

Setting to POLISHING

The standard setting POLISHING is pre-set to WATER OFF and 150RPM.

Settings with a
RotoCom Memory Unit

If RotoPol is connected to RotoCom and communication is in ROTOCOM mode, the following screen will appear:

WATER ON 300RPM
SiC-Paper #320

(Example screen)

- The first line on the screen shows the settings for WATER and SPEED.
- The second line shows the surface to be used, according to the preparation program selected on RotoCom.

DISC  and  toggles between the various surfaces defined in RotoCom.

STRUERS CONSUMABLES

Struers offers a wide range of consumables for both plane and fine grinding. Polishing cloths, diamond and oxide abrasive materials.

RotoPol-11/-15
Instruction Manual

Reference Guide

Table of Contents	Page
1. Advanced Operations	
Selection of Disc	14
Mounting of Specimen Mover (Optional).....	14
Connection to Multidoser (optional)	14
RotoCom Memory Unit (optional).....	14
Setting the Communication	14
Recirculation Cooling Unit (Optional).....	15
Struers Metalog Guide™.....	16
Oxide Polishing.....	16
2. Overview of Menu Screens in the Set-up Menu.....	17
3. Accessories and Connected Equipment	18
4. Trouble-shooting.....	19
5. Maintenance	
Daily Service	20
Weekly Service.....	20
Replacing the Cooling Water.....	20
Checking the Recirculation Cooling Unit.....	20
6. Technical Data	21

1. Advanced Operations

Selection of Disc

RotoPol-11/-15 is designed for MD-Disc of 200 mm dia., wet grinding discs of 230 mm dia. and polishing discs of 200mm dia. For automatic preparation of specimens, only aluminium discs should be used. See the instructions for RotoForce-1.

Mounting of Specimen Mover (Optional)

RotoPol-11/-15 can be equipped with a *RotoForce-1* specimen mover. To mount the specimen mover, see the Instruction Manual of RotoForce-1.

IMPORTANT

If RotoPol-11/-15 has to be relocated, the specimen mover should be removed.

Connection to Multidoser (optional)

RotoPol-11/-15, combined with a specimen mover, can be equipped with *Multidoser* for automatic dosing of suspensions and lubricants. To connect Multidoser, see the appropriate Instruction Manual.

RotoCom Memory Unit (optional)

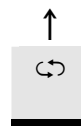
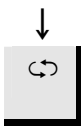
RotoPol-11/-15, combined with RotoForce-1 and Multidoser, can be equipped with a *RotoCom* memory unit. To connect RotoCom to RotoPol, see the appropriate Instruction Manual.

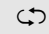
Setting the Communication

- Press MENU  to activate the set-up menu.

COMMUNICATION
STANDARD

Move to the COMMUNICATION menu using the MENU  keys.

Press SELECT  to toggle between STANDARD and ROTOCOM:

COMMUNICATION
ROTCOM

- Select STANDARD if RotoPol is used as a single unit or if combined with a RotoForce-1 and/or Multidoser.
- Select ROTOCOM when RotoPol is connected to a RotoCom unit

- Press STOP  when the desired setting appears.

Recirculation Cooling Unit (Optional)

A Recirculation Cooling Unit (TRECI) can be connected by means of a Recirculation Cooling Unit Connector (ROPRE).

- Connect the cable from the pump in TRECI to ROPRE, according to the diagram of ROPRE.
- Connect the cable with plug from ROPRE to the socket marked Recirculation on the back of RotoPol.
- Connect the cable from ROPRE to the mains.
- Check that the pump rotates in the direction indicated on the pump. If the direction is incorrect, switch two of the phases in the connection to the pump in TRECI (3-phase versions only).
- Replace the outlet tube on the pump in TRECI with the non-return valve with elbow and hexagon nipple from ROPRE. Struers recommends that you seal the threads.
- Connect the pressure hose of RotoPol to the hexagon nipple. Remember to mount the gasket.
- Select the reduction piece from ROPRE which fits the drain hose of RotoPol. Mount the piece on the outlet angle of the TRECI.
- Shorten the drain hose if necessary. Fix the drain hose on the reduction piece: strip away a short length of the steel reinforcement in the drain hose and mount a hose clamp on the stripped end.
- Mount the outlet angle on TRECI.

REMEMBER

Avoid sharp bends in the drain hose and make sure that the hose slopes throughout its length.

IMPORTANT

If a Recirculation Cooling Unit is connected, special precautions must be taken when working with Oxide Polishing Suspensions:

- the water coming from the built-in water tap cannot be used to wash the oxide polishing suspension away. Use clean water.
- The outflow must not be conducted to the Recirculation Cooling Unit. Direct the outlet to the drain or collect the outflow in a separate container.
- When working with Multidoser and Oxide Polishing Suspensions, clean water is used automatically for flushing. Make sure that the outflow is directed to a separate container or to the drain.

Struers Metalog Guide™

RotoPol-11/-15 is designed for preparation of most materials. In Struers Metalog Guide™ you will find a detailed description of grinding/polishing for automated mechanical specimen preparation.

Metalog Guide

Struers Metalog Guide™ offers preparation methods for the most common materials, based on a simple analysis of two key properties: hardness and ductility. Finding the right method is easy, including choice of consumables. Always consult Struers Metalog Guide™ for the correct preparation method for the actual specimens.

Struers Metalog Guide™ contains 6 useful chapters:

Metalogram: a quick and safe guide to the right preparation method.

Metalog Methods: a complete catalogue of preparation methods, based on Struers' vast store of materialographic experience, and employing Struers' range of consumables.

Preparation Philosophy: the basics of modern specimen preparation, seen from a professional point of view.

Metalog Process: the materialographic preparation process from start to finish, logically explained.

Metalog Master: a combined trouble-shooting guide and supply of in-depth information on the processes of mechanical preparation, including an expert system for the solving of preparation problems.

Consumables Specification: quick access to the relevant consumables for the chosen preparation methods.

A complete guide to materialographic specimen preparation.
Contact your local dealer for a free copy of Metalog Guide™.

Oxide Polishing

For polishing with Oxide Polishing Suspension:

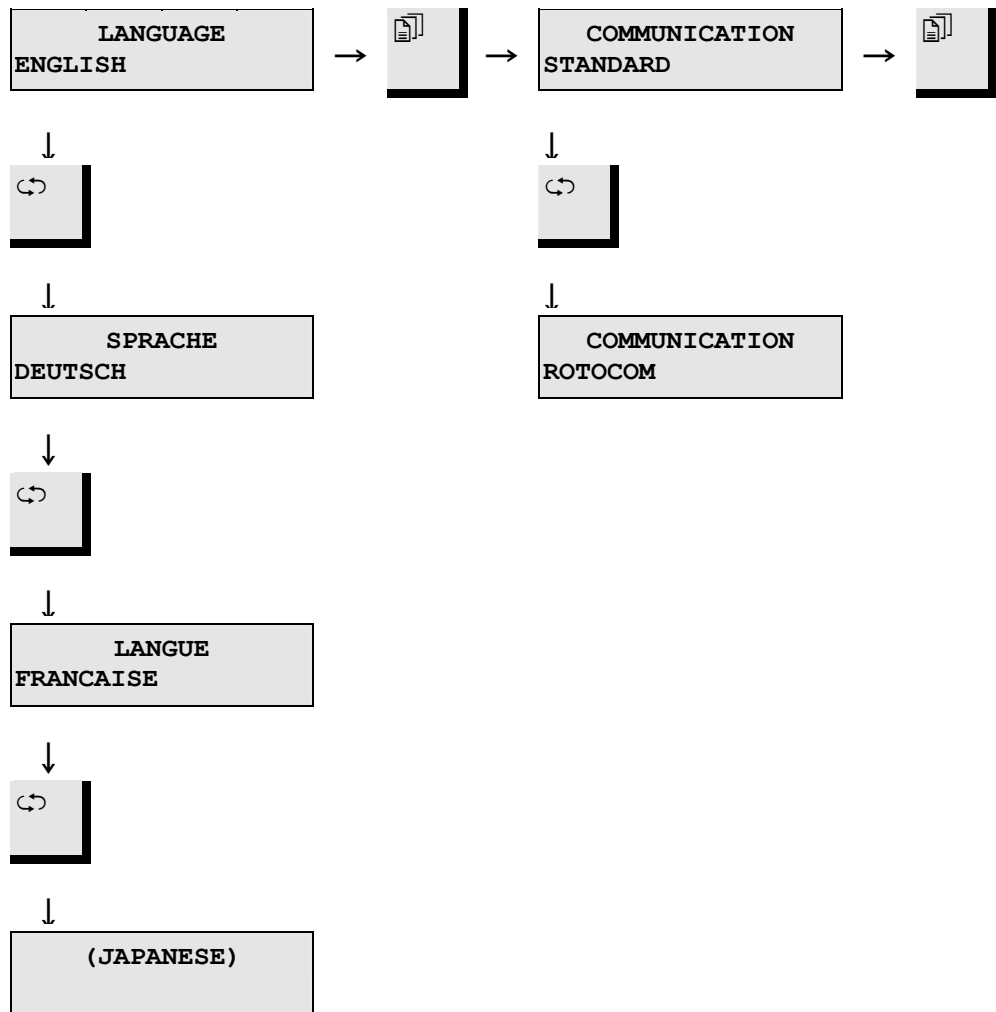
At the end of the polishing process, select WATER ON . Adjust the water tap if necessary. Let the water flow until the oxide polishing suspension has been cleaned away (about 10 seconds).

IMPORTANT

If a Recirculation Cooling Unit is connected, special precautions must be taken:

- the water coming from the built-in water tap cannot be used to wash the oxide polishing suspension away. Use clean water.
- the outflow must not be conducted to the Recirculation Cooling Unit. Direct the outlet to the drain or collect the outflow in a separate container.

2. Overview of Menu Screens in the Set-up Menu


The settings in the top row of the screens show the default values.

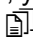3. Accessories and Connected Equipment

Specification	Code
Accessories	
<i>Wet Grinding Discs with Ring</i>	
230 mm dia., PVC	DEPRO
<i>Polishing Discs</i>	
200 mm dia., aluminium	PAMLA
<i>Disc for magnetic fixation</i> 200 mm dia., aluminium	DEMLA
Lid for RotoPol-11/-15	LALID
Connected Equipment	
<i>RotoForce-1, specimen mover</i>	ROFO1
<i>Multidoser (6 Bottles)</i> for automatic dosing of 6 liquids (1-5 suspensions, 0-4 lubricants and 1 OP-Suspension)	MULDO
<i>Multidoser Base Unit (3 Bottles)</i> for automatic dosing of 3 liquids (0-3 suspensions, 0-3 lubricants)	MULBA
<i>RotoCom, memory unit</i>	ROCOM
<i>Recirculation Cooling Unit</i>	TRECI
<i>Recirculation Cooling Unit Connector</i> for connection of Recirculation Cooling Unit (TRECI) to grinding/polishing machines	ROPRE

4. Trouble-shooting

Error	Explanation	Action
Error Message		
OVERLOAD	The motor has been overloaded for a period and the thermo switch has stopped the motor. The motor stops.	Wait until the motor has cooled down. When the display returns to the previous message, the machine can be re-started.
Acoustic Signal		
Long beep.	The selected function is protected. It cannot be activated.	Check the setting.
Machine Problems		
Noise when the machine starts or the machine will not pull.	The belt is not tight enough.	Call a Struers service technician. The belt must be tightened.
Functions are not carried out. RotoCom and/or Multidoser do not operate.	Fuse at the rear of RotoPol blown.	Replace the fuse.
The machine does not operate.	One of the fuses on the right-hand side of the cabinet blown.	Replace the fuse.
Water not draining away.	Drain hose squeezed.	Straighten the hose.
	Drain hose clogged.	Clean the hose.
	Drain hose does not slope downwards.	Adjust the hose to an even slope.
Water dripping underneath the machine.	The water does not drain away.	See above.
Cooling water stops.	Wrong software setting.	Check software setting.
	Water tap on mains closed.	Open for water.
	Built-in water tap closed.	Open for water.
	Built-in water tap blocked.	Clean water tap.
	Filter at the water inlet blocked.	Clean filter.
Cooling water drips after stop.	Defect in the solenoid valve.	Call a Struers service technician. The solenoid valve must be replaced.
The disc vibrates.	Dirt on the underside of the disc or on the turntable.	Clean the contact face between the disc and turntable.
JAPANESE TEXTS	All RotoPol modules have the possibility of showing the screen texts in Japanese. When choosing a European language in the LANGUAGE menu, you might accidentally press  MENU at the Japanese option.	To return to English, German or French texts, just follow this procedure: Press  STOP. Press  MENU. Press  SELECT until you reach the language you prefer. Press  STOP.

5. Maintenance

Daily Service

Clean all accessible surfaces with a moist cloth.

Weekly Service

Remove the grinding/polishing disc and remove all dirt from the tray and the drain tube.

Replacing the Cooling Water

Only relevant when a Recirculation Cooling Unit is connected: replace the cooling water in the Recirculation Cooling Unit at least once a month.

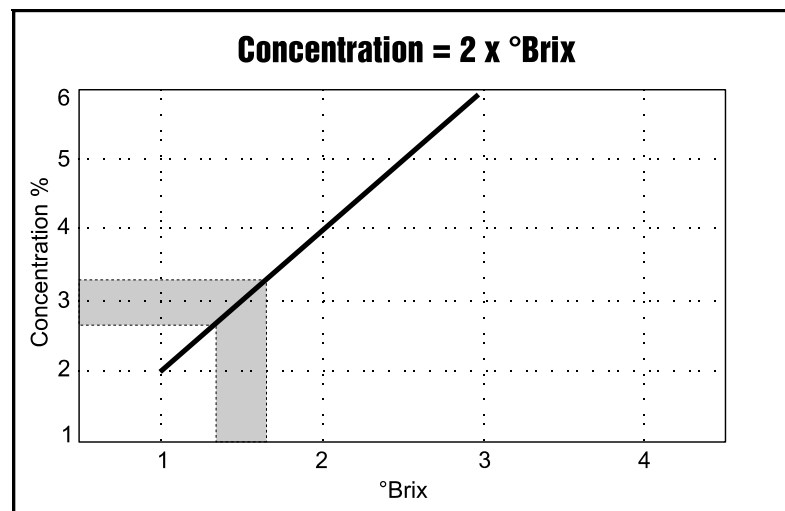
- Place a Disposable Plastic Container (TREPO) in the tank and fold it over the edge.
- Fill the tank with 30 l of water. NB! Too high a level of coolant in the tank could damage the pump.
- Add Struers Additive according to the instructions.
- Put the lid on the tank.
- Place the pump in the lid

IMPORTANT

Always maintain the correct concentration of Struers Additive in the cooling water (percentage stated on the container of the Additive). Remember to add Struers Additive each time you refill with water.

Checking the Recirculation Cooling Unit

The cooling unit should be checked for cooling water every week and refilled when the flushing pump cannot reach the cooling water. Remember to add Struers Additive for Cooling Fluid: One part of Additive for 33 parts of water. The concentration of additive should always lie between 2.7 and 3.3%; the pH should be around 9. Replace the cooling water if pH is ≤ 8.0 . To check the concentration of additive, use a refractometer (0-10 Brix, see diagram.). Concentration = $2 \times \text{Brix}$. Add Struers Additive for Cooling Fluid if the concentration is lower than 2.7%. Add water if the concentration is higher than 3.3%.

6. Technical Data

Subject		Specifications		
		RotoPol-11	RotoPol-15	
Disc	Diameter	200-230 mm / 8-9"		
	Speed	150/300 rpm	40-600 rpm	
	Rotational direction	Anti-clockwise		
	No. of motors	1		
	Motor	90W / 0.12HP at 150 rpm		
		180W / 0.24HP at 300 rpm	370W / 0.49HP at 600 rpm	
	Torque at disc	Continuous 8Nm / 5.9lb-ft	Continuous 6Nm / 4.4lb-ft at 600 rpm	
Max. 13Nm / 19.6lb-ft		Max. 12Nm / 18.1lb-ft		
Connected specimen movers	RotoForce-1	1		
Automatic stop	RotoForce-1 and Multidoser	Yes		
Data exchange	RotoCom	Yes		
Software and electronics	LC display	2x16 characters		
	Controls	Touch pads		
	Memory	EPROM/RAM/NV-RAM		
Fittings	No. of water taps	1		
	Splash rings	1		
Safety standards	General	89/392/EEC & 91/386/EEC & 93/44/EEC EN292-1, EN292-2, EN60204-1 (IEC204-1)		
	EMC	89/336/EEC & 92/231/EEC EN50081-1, EN50082-1 FCC part 15, subpart B		
Noise level	At idle running, at a distance 1.0 m / 39.4" from the machine	52 dB (A)	55 dB (A)	
Surrounding temperature		5-40°C / 41-104°F		
Humidity	Non condensing	0-95%RH		
Supply	Power	1 phase (N+L+ PE)		
	Power consumption	360W / 0.48HP	580W / 0.60HP	
	Voltage / frequency	Max continuous load		
	1 x 100V / 50-60Hz	6.0A	6.0A	
	1 x 110-120V / 50-60Hz	6.0A	6.0A	
	1 x 220-240V / 50-60Hz	3.0A	3.0A	
	Pressure for tap water	1-10 bar / 14.5-145 psi		
	Water inlet	1" or 3/4"		
	Water outlet	ø32 mm / 1 1/4"		
Dimensions and weight	Width	410mm / 16.1"		
	Depth	670mm / 26.4"		
	Height	315mm / 12.4"		
	Weight	22Kg / 48.5lbs		

Quick Reference Guide

Mounting a Disc

Place the disc on the turntable and turn it until the three driving pins engage with the holes in the turntable.

Starting the Disc

Press START . The disc starts rotating. The selected functions are executed and connected RotoSystem modules start operating.

Stopping the Disc

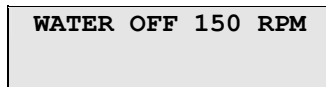
Press STOP . The disc stops rotating. The selected functions and connected RotoSystem modules are stopped.

Grinding/Polishing


- Mount the required surface on a disc.
- Mount the disc on the turntable.
- Select the desired functions, if RotoCom is not connected.
- Press START .


Settings

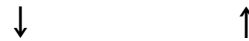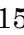Setting to User Select

The User Select screen has an empty line at the bottom.

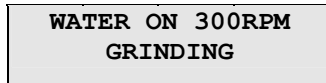
- Press  or  to select WATER ON or WATER OFF.
- Press SPEED  or  to select the speed. (RotoPol-11)
- Press SPEED  or  to adjust the speed. (RotoPol-15)

Press DISC  or  to change between the settings.



Setting to GRINDING

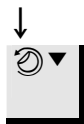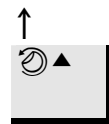
The standard setting GRINDING is pre-set to WATER ON and 300RPM.

Press DISC  or  to change between the settings.

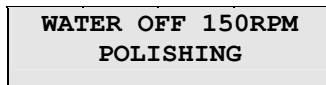Setting to POLISHING

The standard setting POLISHING is pre-set to WATER OFF and 150RPM.

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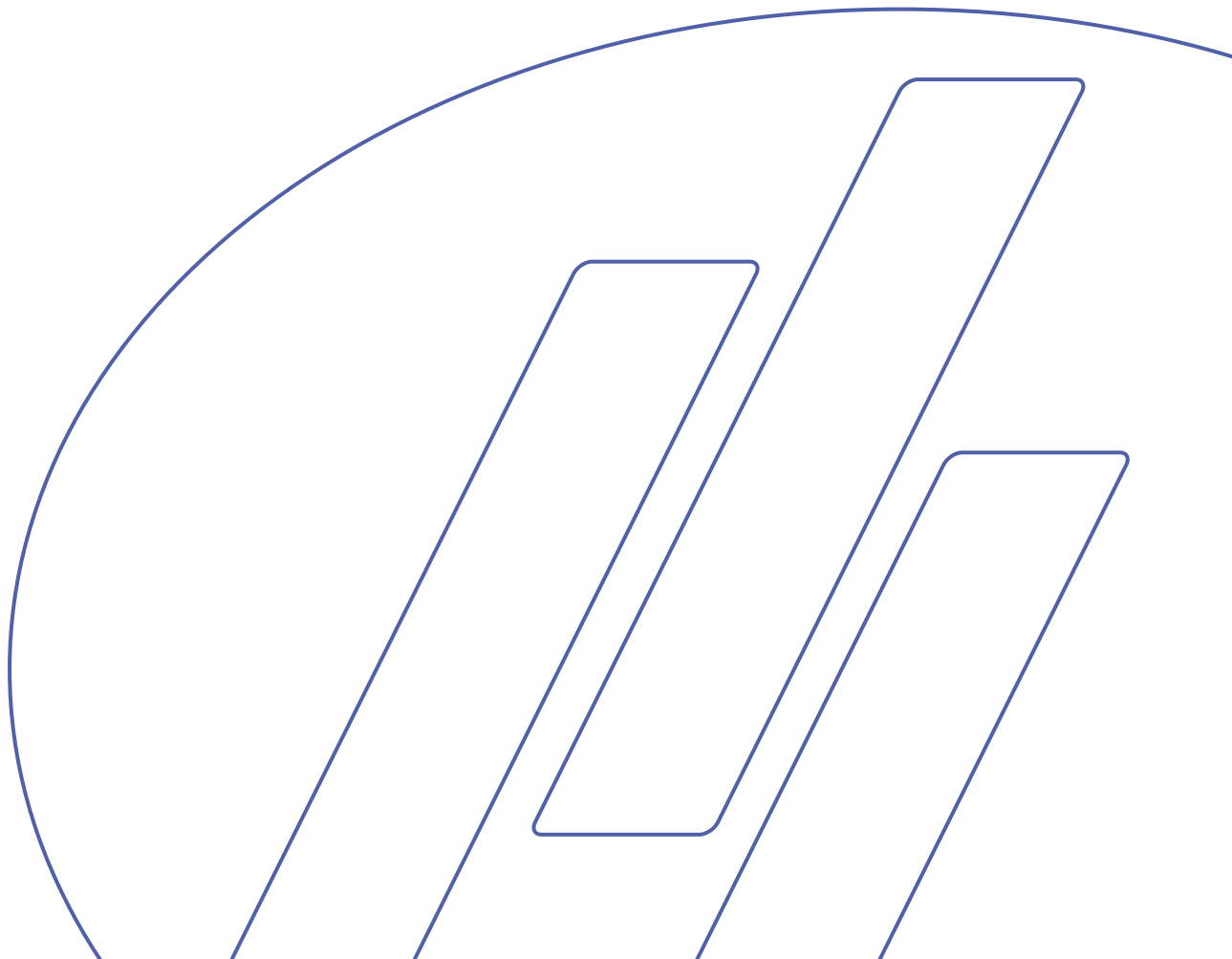
Spare Parts and Diagrams

RotoPol-11/-15
Spare Parts and Diagrams

**Always state *Serial No* and *Voltage/frequency*
if you have technical questions or when ordering spare parts.**

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
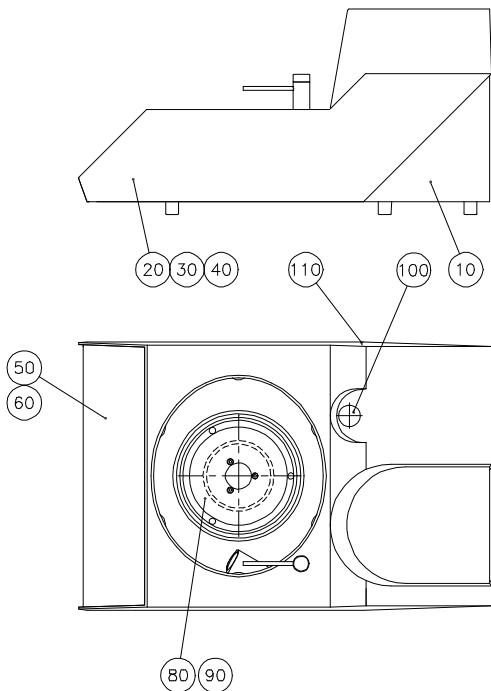
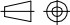
Some of the drawings may contain position numbers
not used in connection with this manual.

RotoPol-11/-15
Spare Parts and Diagrams


Drawing	Pos.	Spare Part	Part No.
15190002	100	Cover plate	2GT62740
	110	Cover plate	2GT61309
14910061		<i>Water Tap, Complete</i>	14910061
		Gaskets for water tap	14910071
	090	Water tube and tap, complete	14910062
<i>Various Parts</i>		Pressure hose, white	359MP001
		Connection piece 1/2"	2NG30013
		Gasket Ø11/Ø24 x 1.5, (10 pcs.)	R3590359
		Gasket 3/4" with filter, (10 pcs.)	445MP068
		Danflex K-126 Ø32, incl. worm belt	RNU30232
		Protection ring Ø250	15180101
		Insex key 5	2GR00050
RotoPol-11 only			
15190022	40	Poly-V belt, (Length 1016 mm)	2JD31016
	270	Fuse 10.0A (100-120V), (10 pcs.)	372MP075
	270	Fuse, 6.3A (220-240V), (10 pcs.)	315MP134
	310	Fuse, 2.5A, (10 pcs.)	RFU14000
RotoPol-15 only			
15260010	40	Poly-V belt, (Length 965 mm)	2JD30965
	270	Fuse 10.0A, (10 pcs.)	372MP075
	310	Fuse, 3.15A, (10 pcs.)	392MP076

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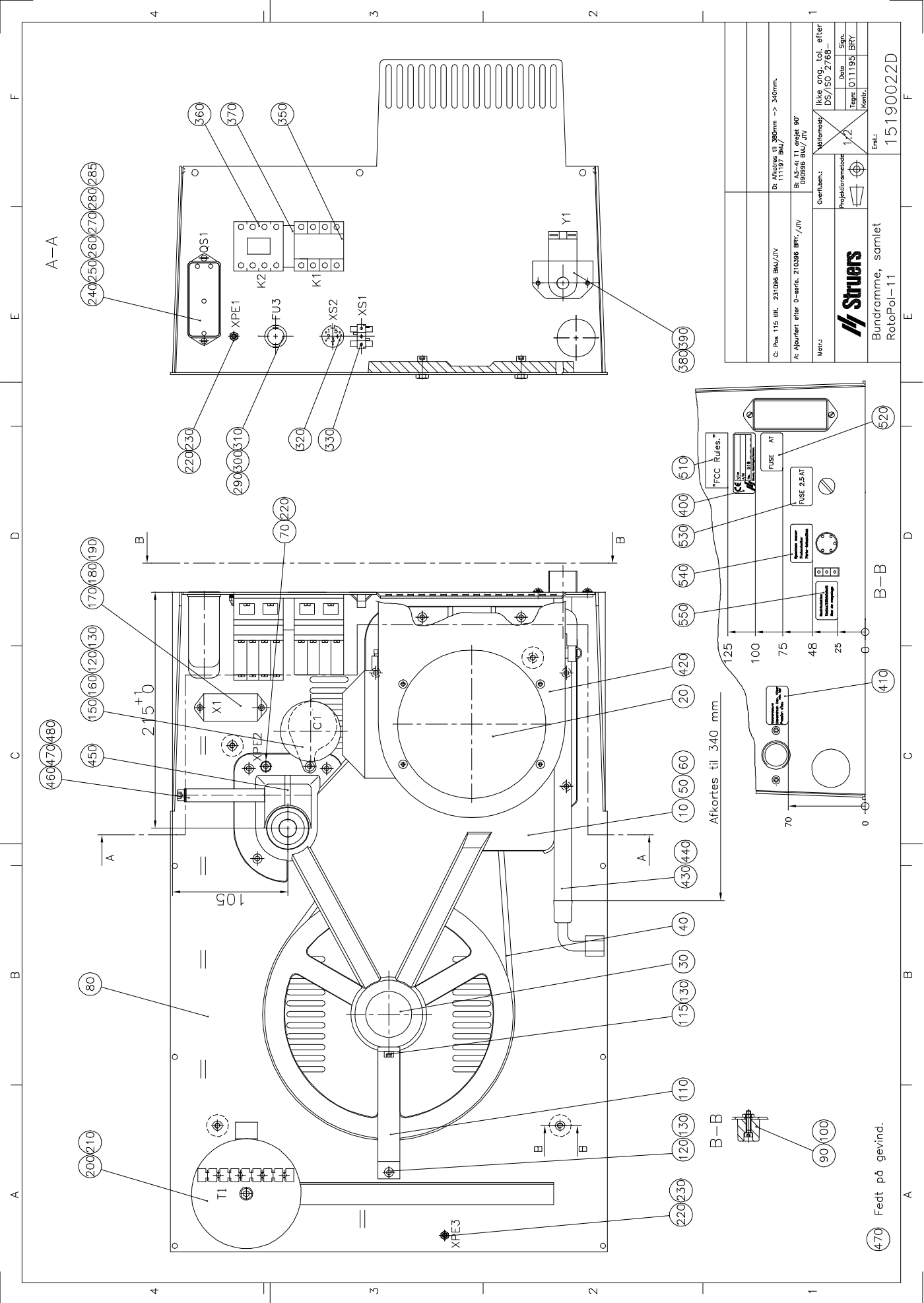
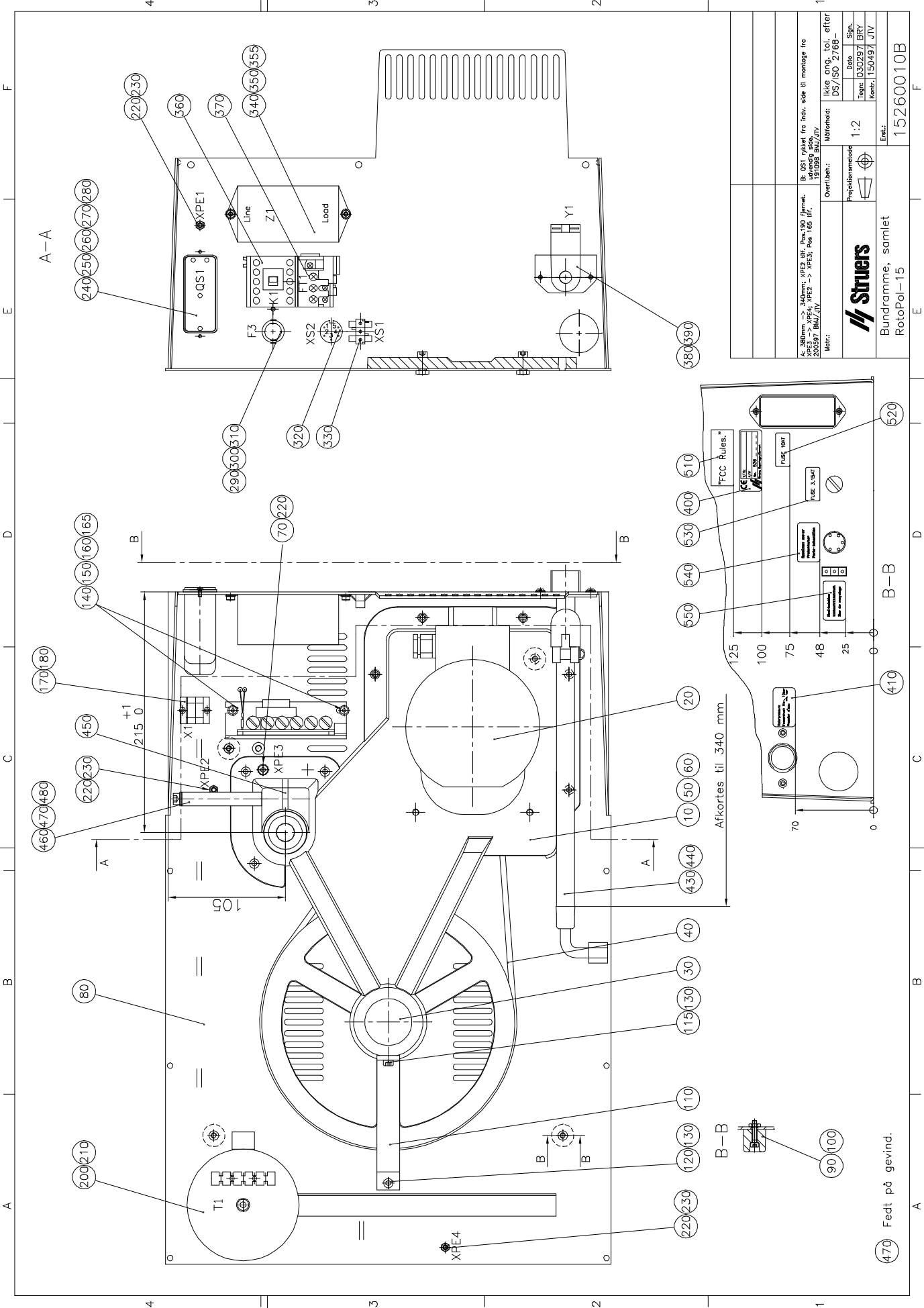
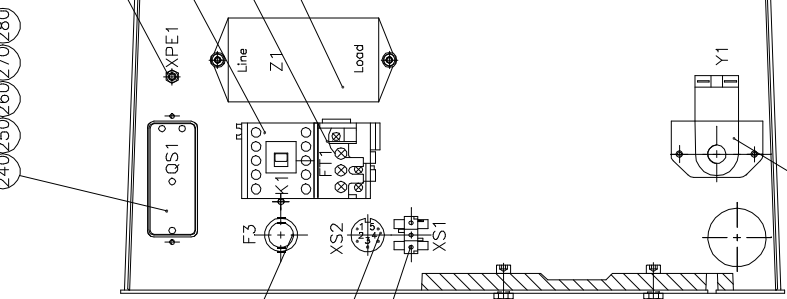
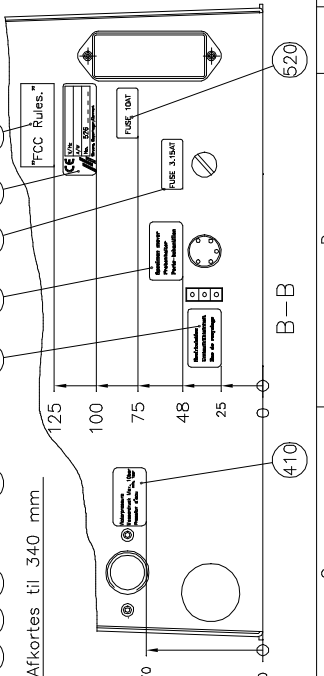
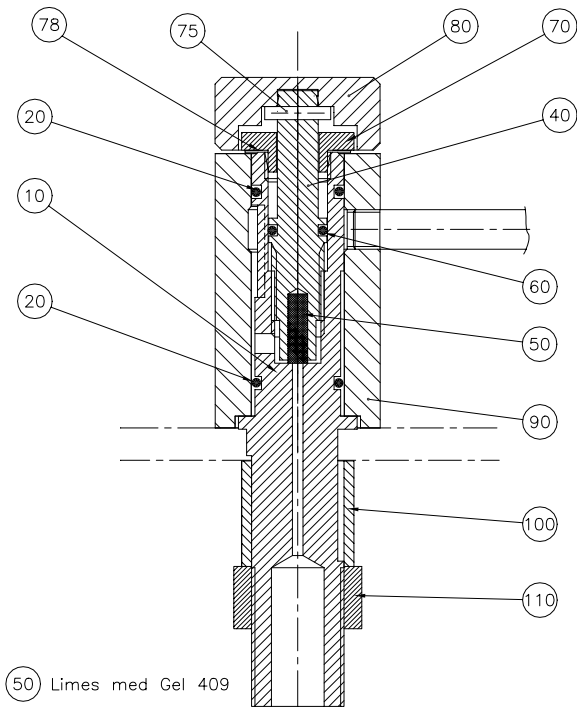
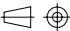
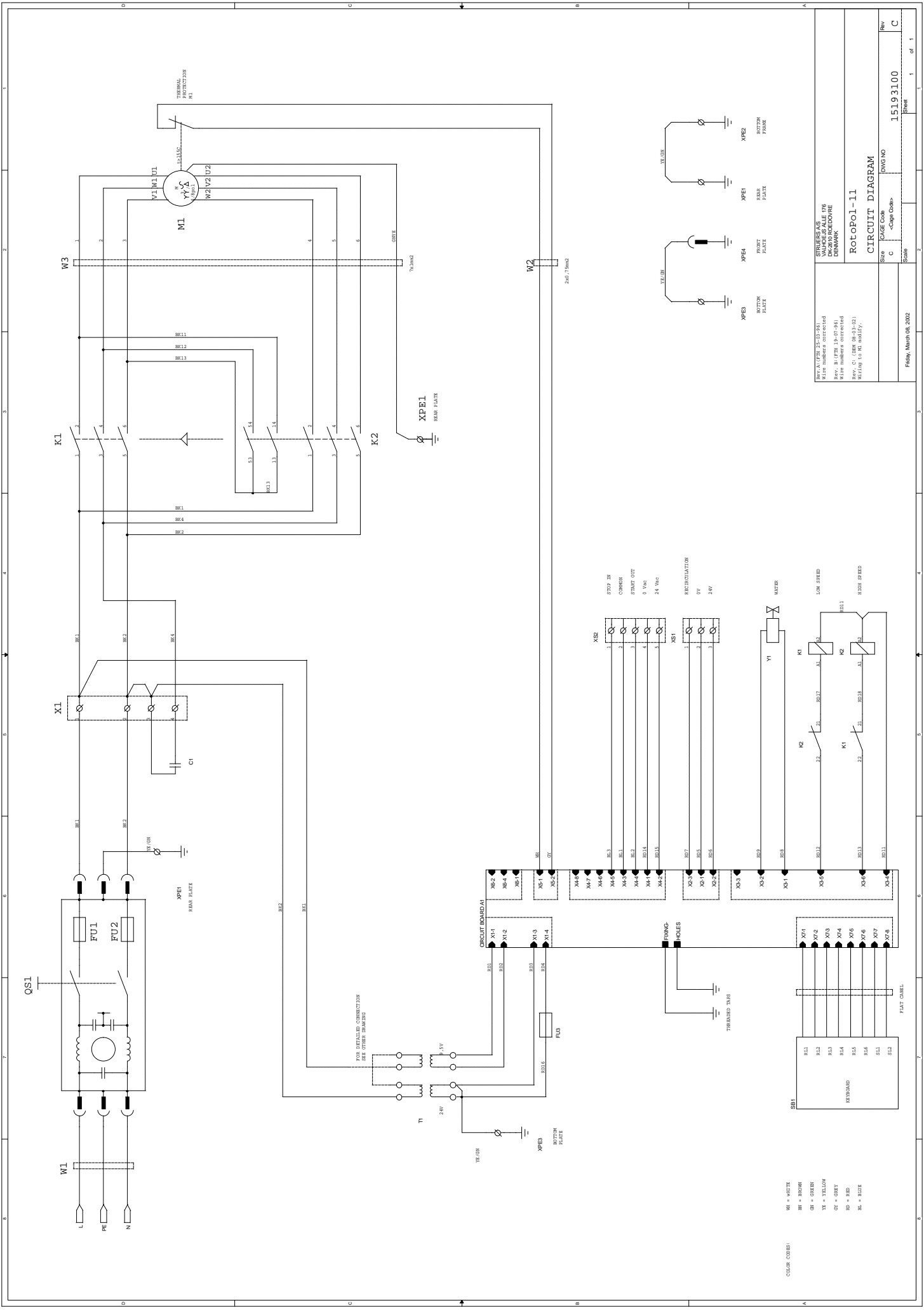
Scale: 1:1
 DWG NO: 15193100
 Rev: C

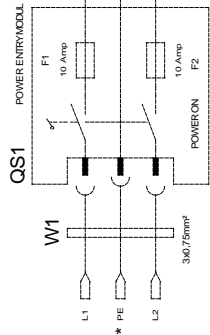Friday, March 08, 2002


* MUST BE SHORTCUT
FIRST COIL WINDING
REGULATIONS
F = 15 Amp MAXIMUM

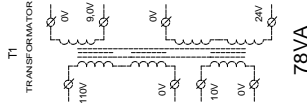FOR DETAILED CONNECTION
SEE DRAWING 1572119












78VA

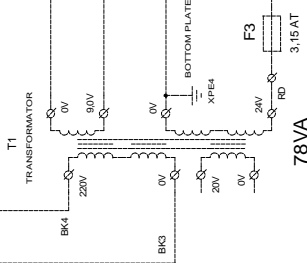FOR DETAILED CONNECTION
SEE DRAWING 1492319

78VA

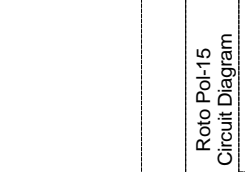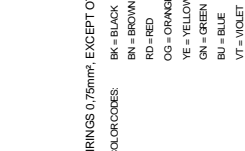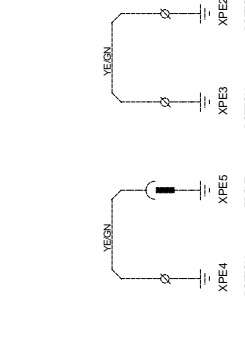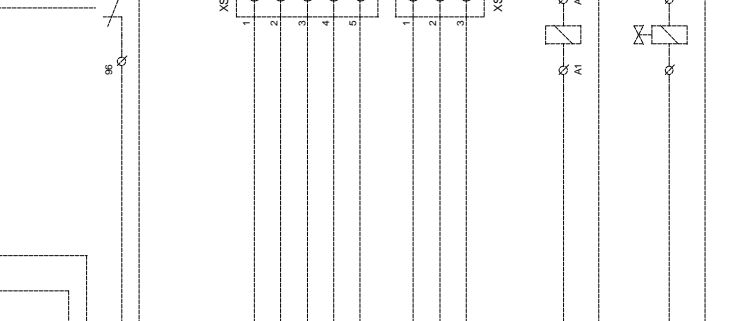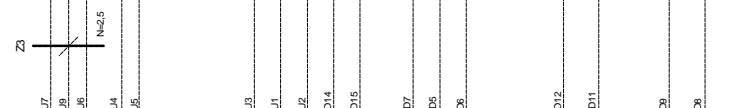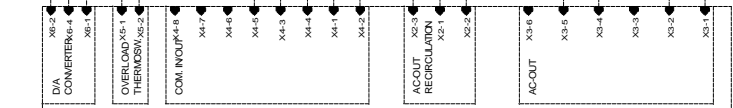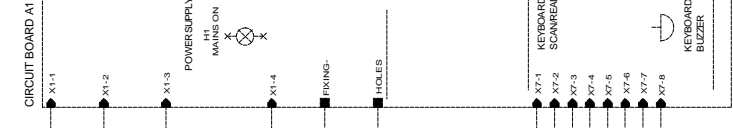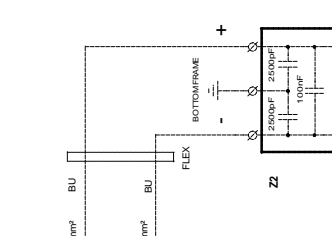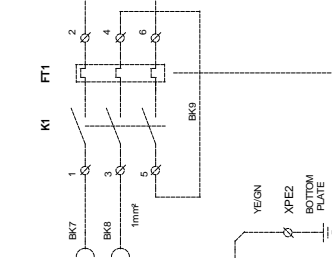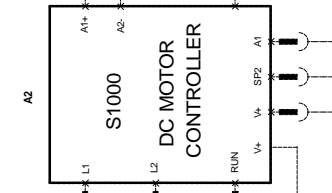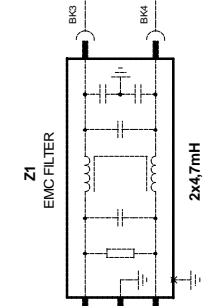ALL WIRINGS 0.75mm², EXCEPT OTHERWISE MARKED.


- COLORCODES:
- BK = BLACK
 - BN = BROWN
 - RD = RED
 - OG = ORANGE
 - YE = YELLOW
 - GN = GREEN
 - BU = BLUE
 - VT = VIOLET
 - GY = GREY
 - WH = WHITE

STRUERS A/S KVLAVSVEJ 10 DK-2600 SØNDERBY DENMARK TELEPHONE: 33203500	Size A2	CAGE CODE SCALE	DWG NO 15263100	Sheet 1 of 1
---	------------	--------------------	--------------------	-----------------

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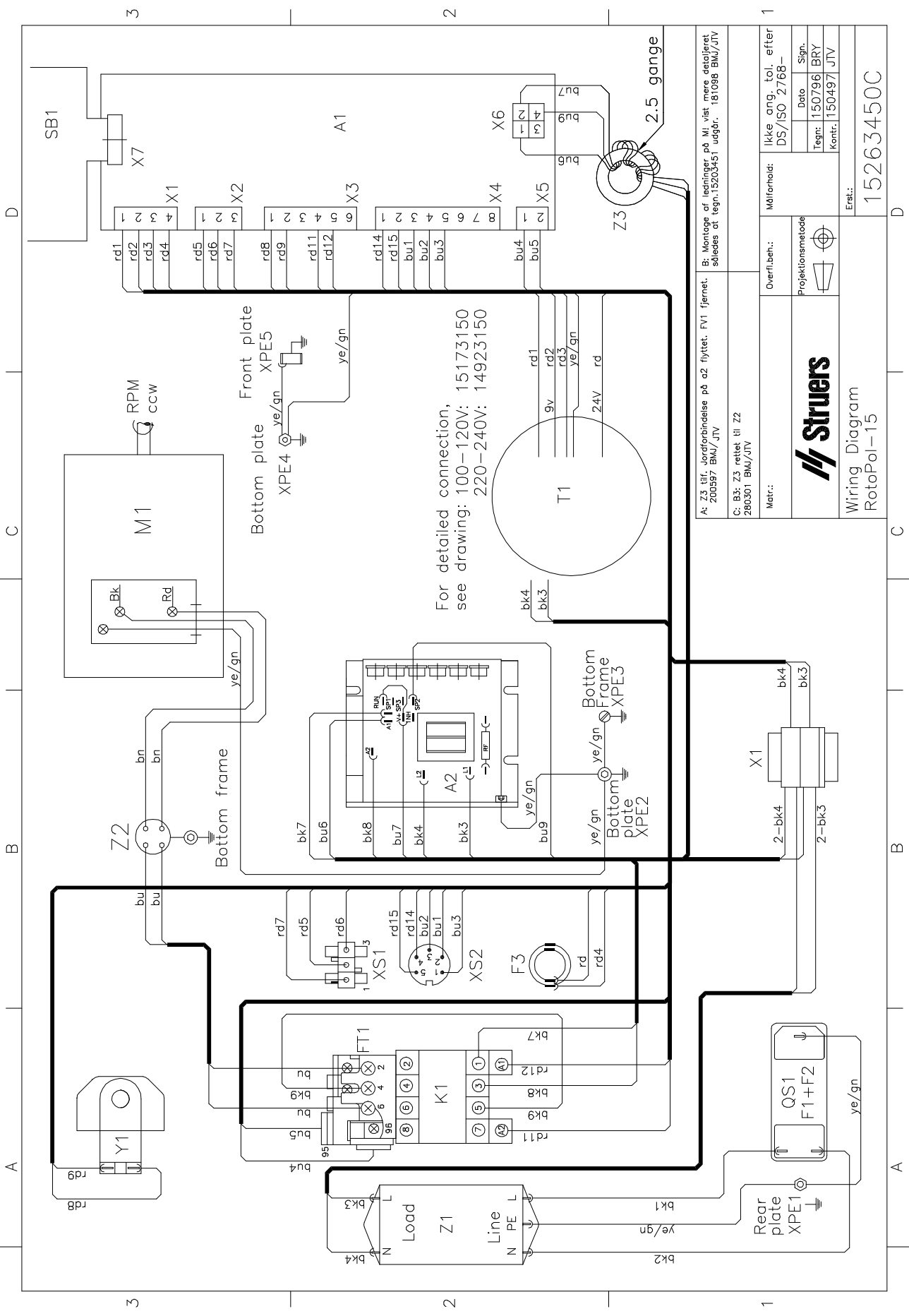Herewith declares that

<i>Product Name:</i>	RotoPol-11, -15
<i>Type No:</i>	519, 526
<i>Machine Type:</i>	Grinding and polishing machine

is in conformity with the provisions of the following directives:

Safety of Machinery 89/392/EEC and 91/368/EEC and 93/44/EEC according to the following standard(s):
EN292-1:1991, EN292-2:1991, EN60204-1:1992.

EMC-Directive 89/336/EEC and 92/31/EEC according to the following standard(s):
EN50081-1:1991, EN50082-1:1992.

Low Voltage Directive 73/23/EEC and 93/68/EEC according to the following standard(s):
EN60204-1:1992.

Supplementary Information The equipment complies with the American standards:
FCC part 15, subpart B and UL508.

The above has been declared according to the global method, module A

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Klaus Kisbøll, Entwicklungsleiter, Struers A/S
Klaus Kisbøll, Chef du service développement, Struers A/S